QUANTOFIX® Arsen 10

de

Inhalt:

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en

Pack content:

- | | |
|--|-------------------|
| 1 aluminium container with 100 test strips | 1 reaction vessel |
| 1 bottle of As-1 | 1 syringe 10 mL |
| 2 bottles of As-2 | 1 needle |
| 2 bottles of As-3 | |

Safety precautions:

Each test field on the test strips contains mercury(II)bromide. Toxic in contact with skin. As-2 contains sulfamic acid 50-100 % CAS 5329-14-6. WARNING.

General indications:

Remove only as many test strips as are required. Close the container immediately after removing a strip. Do not touch the test field.

Instructions for use:

- Using the syringe, fill 3 x 10 mL sample solution into the reaction vessel. After adding each of the following reagents, shake gently for about 2 s:
- 1 drop of As-1
- 1 measuring spoon of As-2
- 1 measuring spoon of As-3
- Immediately afterwards dip the test strip with the test field about 2 cm deep into the reaction vessel and then fit the lid (see drawing). To vent the reaction vessel, pierce the lid with the needle and leave the needle in the lid.
- Wait 10 minutes.
- Afterwards dip the test strip with the test field into distilled water for two seconds.
- Shake off excess liquid and compare with the color scale.

Interferences:

The following ions interfere with the determination only when the concentrations below are exceeded:

1000 mg/L: Ca²⁺, K⁺, Mg²⁺, Na⁺, Zn²⁺, Cl⁻, CN⁻, CO₃²⁻, SO₄²⁻, EDTA;
 500 mg/L: Fe²⁺, Fe³⁺, F⁻; 100 mg/L: Al³⁺, Sn²⁺, CrO₄²⁻, NO₂⁻, NO₃⁻, PO₄³⁻; 25 mg/L: ClO₃⁻;
 5 mg/L: Co²⁺; 2 mg/L: S²⁻, SO₃²⁻; 1 mg/L: Ni²⁺, Sb³⁺, SeO₃²⁻, S₂O₃²⁻; 0.5 mg/L: Ag⁺, Cu²⁺

Disposal:

Dispose of used strips at an authorised hazardous waste collection point.

Storage:

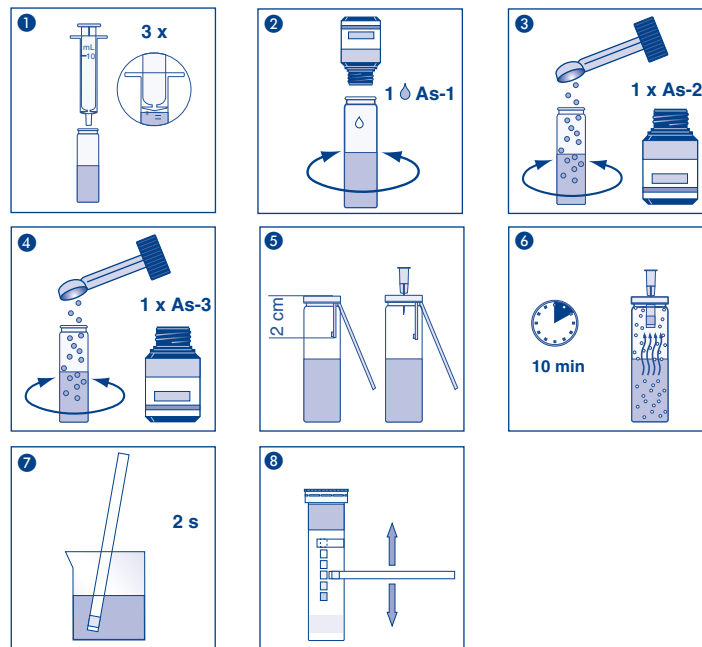
Avoid exposing the strips to sunlight and moisture. Keep container cool and dry (storage temperature 4–30 °C)..

If correctly stored, the test strips may be used until the use-by-date printed on the packaging.

Additional information:

The test strip container stopper contains a non-toxic drying agent. If swallowed, drink plenty of water.

MACHEREY-NAGEL GmbH & Co. KG · Valenciener Str. 11 · 52355 Düren · Germany
 Tel.: +49 24 21 969-0 · info@mn-net.com · www.mn-net.com

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MACHEREY-NAGEL GmbH & Co. KG · Valenciennr Str. 11 · 52355 Düren · Germania